



Day 1: Immunohistochemistry – TDP-43 Aptamer

Things to do while you're waiting for your timer

- ✓ Get the blocking reagents from the cold room

- ✓ Make up 10mM citric acid retrieval agent 1:10 from stock.
Stock: 100 mM citric acid (19.12g in 1L dH₂O).
- ✓ Make up 1:10 TBS (Tris:saline)
- ✓ If using, dilute citric acid retrieval 1:10 (stock: dH₂O)

The steps

Deparaffinise slides

Xylene	3 min	<input type="checkbox"/>
Xylene	3 min	<input type="checkbox"/>
Alcohol	2 min	<input type="checkbox"/>
Alcohol	2 min	<input type="checkbox"/>

Place slides in tap water (but see below).

NB: Do not place slides in tap water if employing optional picric acid step below

(Optional step) Remove Formalin Pigment

Saturated Alcoholic Picric acid	15 min	<input type="checkbox"/>
Wash in warm, running tap water	15 min	<input type="checkbox"/>

Antigen retrieval

Put 500ml dH ₂ O in the pressure cooker before starting	15 min	<input type="checkbox"/>
--	--------	--------------------------

Coverplate

dH ₂ O	5 min	<input type="checkbox"/>
Peroxidase block (3 drops)	30 min	<input type="checkbox"/>
Wash with TBS	5 min	<input type="checkbox"/>
Avidin block (4 drops)	15 min	<input type="checkbox"/>
Wash with TBS	5 min	<input type="checkbox"/>
Biotin Block (4 drops)	15 min	<input type="checkbox"/>
Wash with TBS	5 min	<input type="checkbox"/>
Make up aptamer in MilliQ H ₂ O		<input type="checkbox"/>
Wash with dH ₂ O	5 min	<input type="checkbox"/>
Aptamer incubation (1in500)	3 hours	<input type="checkbox"/>
Fix step 4% PFA in dH ₂ O	Overnight in fridge	<input type="checkbox"/>



Day 2: Immunohistochemistry – TDP-43 Aptamer

	Re-coverplate (optional)		
	Wash with dH ₂ O	5 min	<input type="checkbox"/>
✓ Make up DAB Chromogen (1:20)	Anti-Biotin/HRP 1:100 in MilliQ	30 min	<input type="checkbox"/>
1 µl DAB chromogen + 19 µl DAB substrate buffer	Wash with dH ₂ O	5 min	<input type="checkbox"/>
So ...			
50 µl DAB chromogen + 950 µl DAB substrate buffer	DAB (1:20)	5 min	<input type="checkbox"/>
	Wash with dH ₂ O	5 min	<input type="checkbox"/>
	Wash well in running tap water		<input type="checkbox"/>
	Counterstain with haematoxylin		<input type="checkbox"/>
	Blue in lithium carbonate		<input type="checkbox"/>
	Dehydrate, clear and mount		<input type="checkbox"/>

References for citation of this method

Please cite both of these if using this method:

The citation, Spence and Waldron *et al.*, 2024, for the first publication for the development, modification and employment of the TDP-43 RNA aptamer to stain human tissue published in *Acta Neuropathologica*.

RNA aptamer reveals nuclear TDP-43 pathology is an early aggregation event that coincides with STMN-2 cryptic splicing and precedes clinical manifestation in ALS.

Holly Spence*, Fergal M. Waldron*, Rebecca S. Saleeb, Anna-Leigh Brown, Olivia M. Rifai, Martina Gilodi, Fiona Read, Kristine Roberts, Gillian Milne, Debbie Wilkinson, Judi O'Shaughnessy, Annalisa Pastore, Pietro Fratta, Neil Shneider, Gian Gaetano Tartaglia, Elsa Zacco, Mathew H. Horrocks, Jenna M. Gregory[†]. 2024. *Acta Neuropathologica* 2024 Mar 5;147(1):50. DOI: 10.1007/s00401-024-02705-1. *equal contributions, [†]corresponding author.

The citation for the SOP and tick-sheet, Waldron and Spence *et al.*, 2024, published on *protocols.io* is:

TDP-43 RNA aptamer staining to detect pathological TDP-43 in FFPE human tissue, as described in Spence and Waldron *et al.*, 2024 (*Acta Neuropathologica*): A SOP and tick-sheet. v2.

Fergal M. Waldron*, Holly Spence*, Jenna M. Gregory[†]. *protocols.io* 2024; DOI: dx.doi.org/10.17504/protocols.io.eq2lyjo4mlx9/v2. *equal contributions, [†]corresponding author.