

eDNA[®] citizen SCIENTIST

ASSEMBLY AND USE INSTRUCTIONS




The Citizen Scientist uses Smith-Root self-preserving filters:

Part #	Description
11579	Filter Pack, 1.2µm Self-Dessicating
11580	Filter Pack, 5µm Self-Dessicating
11747	Filter Pack, .45µm Self-Dessicating

Items manufactured by companies other than Smith-Root carry the original manufacturer's warranty. Please contact product manufacturer for return instructions.


All Smith-Root, Inc. manufactured products are covered by a one year warranty.

Credit & Refund Policy: Customers returning equipment, in new condition, will be given credit five days from the date of the return. A return authorization must accompany returns. Valid equipment returns include, but are not limited to, ordering incorrect equipment, funding deficits, and defective equipment returned for reimbursement. All returns are subject to a restocking fee and applicable shipping charges. The restocking fee is figured at 10% of the purchase price but not less than \$20.00. Customers receiving equipment in damaged condition will be referred to the shipping company for insurance reimbursement.

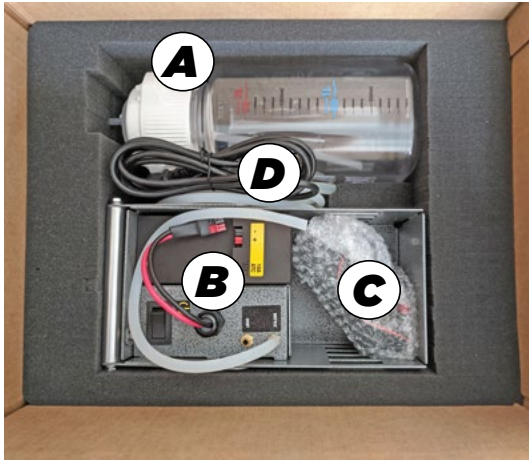


PACKAGE CONTENTS

- A. Vacuum bottle
- B. Caddy with pump and battery
- C. Battery charger
- D. Battery charger power cable

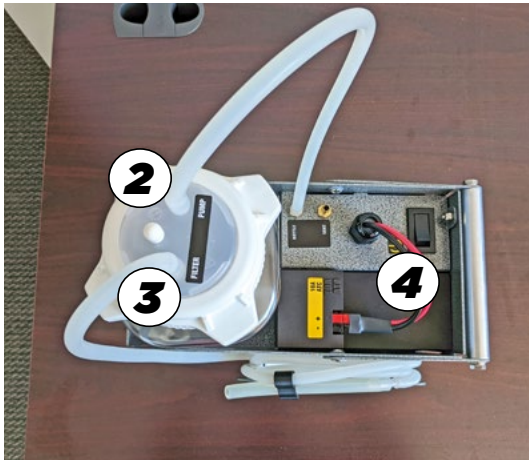

ASSEMBLY INSTRUCTIONS

1. Place bottle in caddy
2. Connect pump to bottle
3. Connect bottle to filter tube
4. Connect power cable
5. Elevate caddy handle

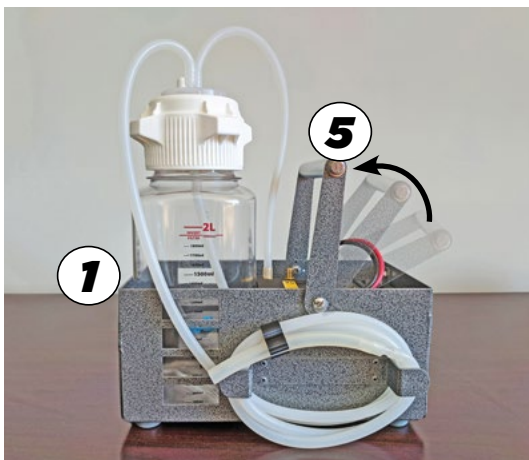

You are ready to sample!

INSTRUCTIONS



1

Find a level spot



2

Open filter packet



3

Attach snorkel tube



4

Connect bottle to filter



5

Turn on vacuum pump



6

Insert snorkel in water



Invert filter at water line



Air dry filter 20 seconds



Turn off vacuum pump



Place filter in zip-top bag



Seal bag shut and label



Discard filtrate water



info@smith-root.com
(360) 573-0202
Vancouver, WA USA
www.smith-root.com

